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C.K. YEUNG WORLDWIDE LIMITED

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We delivers customer growth & productivity ...

\$1.7B Revenue

10,700 employees

Operating in 125 countries

Member S&P 500 Index

Division : Life & Analytical Science

Division: Opto-Electronics

Global technology leader providing products and services to customers in health sciences and other advanced technology markets that require innovation, precision and reliability

FROST & SULLIVAN

PerkinElmer, Inc.

Analytical Instrumentation Company of the Year

2003

PerkinElmer AcycloPrime-FP SNP Detection System Named 'Product of the Year' by Genome Technology Magazine

BOSTON, Oct 7, 2002 (BUSINESS WIRE) -- PerkinElmer, Inc. (NYSE: PKI), a leading provider of drug discovery, life science research and genetic disease screening solutions, was selected for "Product of the Year" honors for its AcycloPrime(TM)-FP SNP Detection System by readers of Genome Technology magazine, an independent provider of news and information covering the business and technology of genomics, proteomics, and bioinformatics. PerkinElmer was also a finalist in the "Company of the Year" category. The Company was honored and presented a trophy at a ceremony recognizing the 2002 Genome Technology Leaders at The Institute for Genomics and Bioinformatics, held in Boston.

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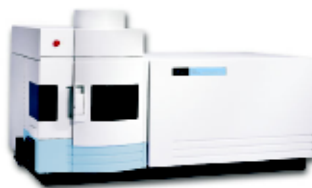
Email through@innovation,precisionandreliability.com

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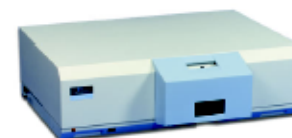
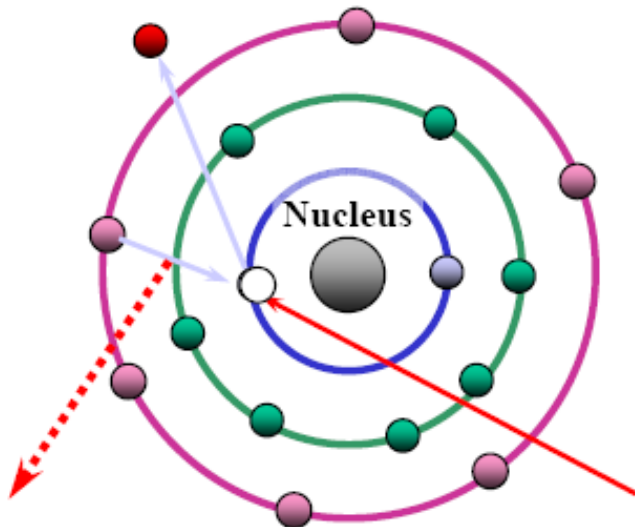
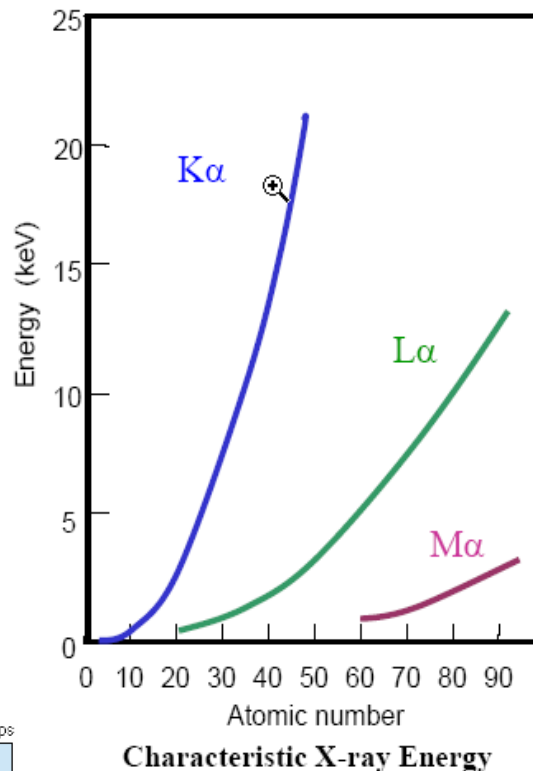
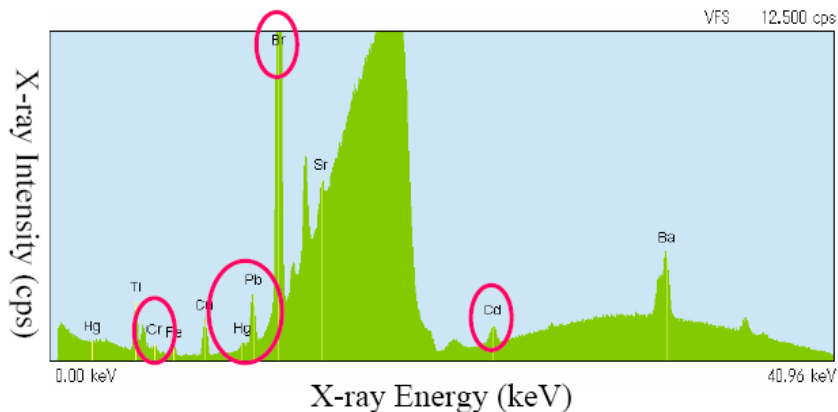
Fluorescent X-ray

Primary X-ray

X-ray Energy (keV) → Element
 X-ray Intensity (cps) → Concentration

Ltd. Tel: 2802-3322 Fax: 2802 8988

Email: joseph@ckyeung.com

Why EDXRF is a popular tool in metal analysis ?

- Fast (real time information display), easy to operation
- Excellent screening tool for metals; 1.0 – xx%
- Non-destructive sample pre-treatment
- Low maintenance

XRF 之分析方法 *Disadvantage*

➤ Measurement time 1,800 sec.

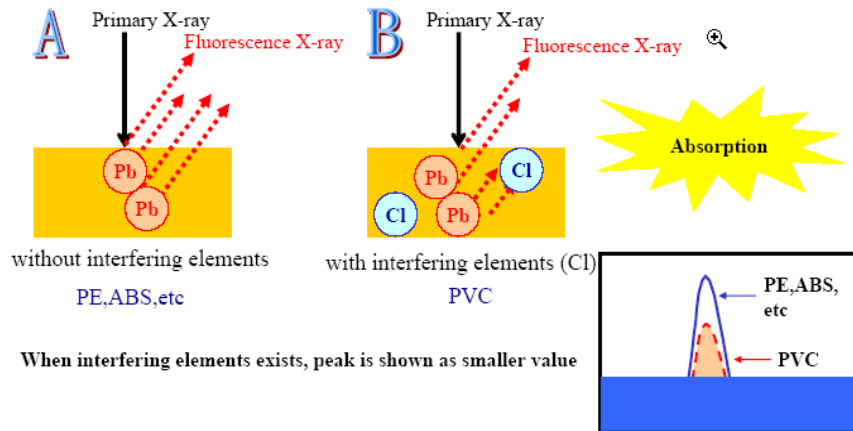
- Detection limit Pb, Cd, Br and Hg : 2 ppb
- Cr : 5 ppm

➤ Measurement time 100 sec.

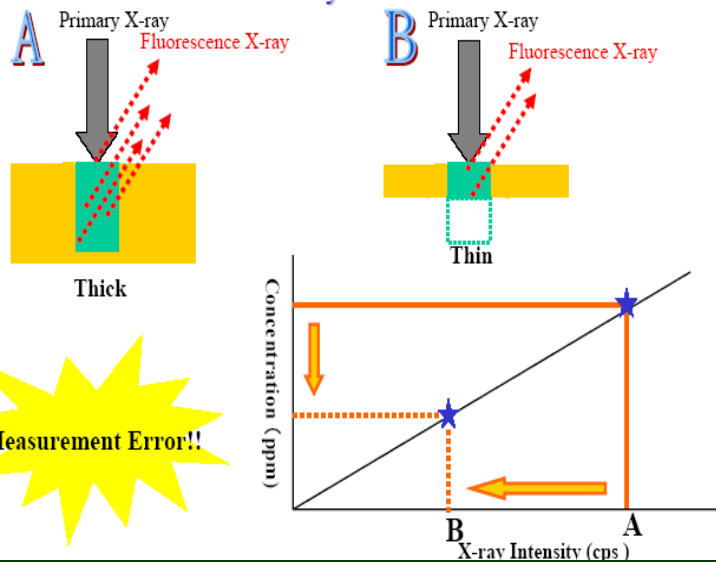
- Detection limit Pb : 36 ppm
- Cd : 20 ppm
- Br and Hg : 40 ppm
- Cr : Approx. 100 ppm

Sample: PVC

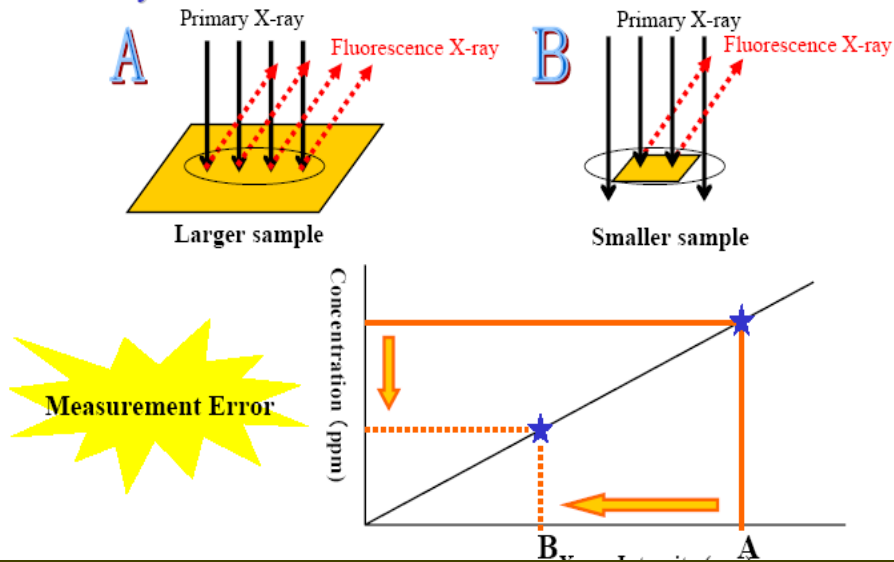
■ Absorption of Cl

➤ Excitation volume of X-rays

➤ X-ray irradiation area

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SECURITY GRADE
CLASS B

MANUAL FOR MEASUREMENT OF Cd AND Pb
BY NITON XRF (XL-700 SERIES) ANALYZER

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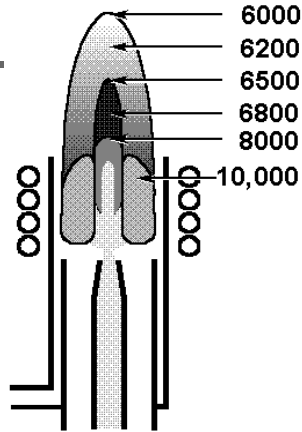
Temperature (K) $\pm 10\%$

Absorption (ν_a)

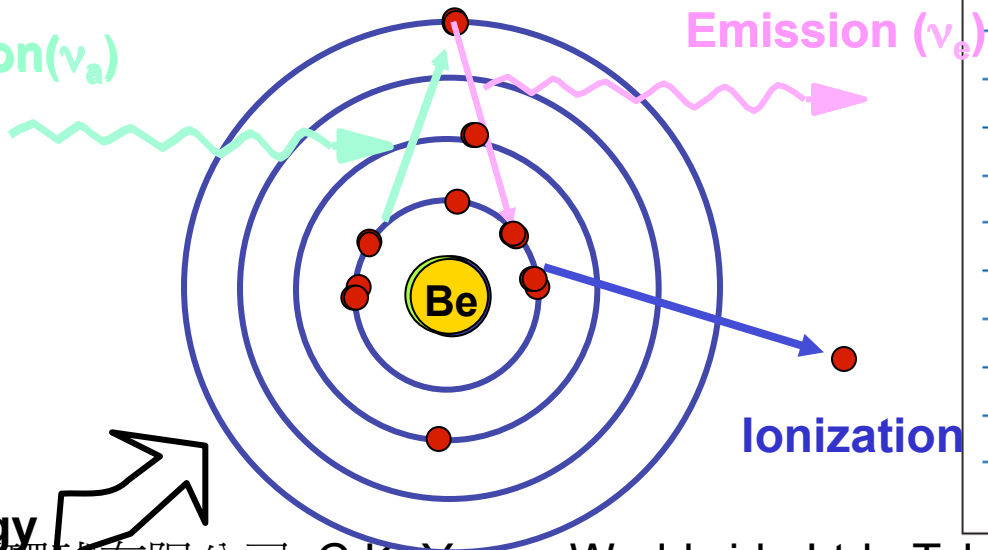
Table 1. Typical Optima 2100 DV detection limits in $\mu\text{g/L}$ (ppb) based on three standard deviations. All values were obtained using axial viewing, a GemCone nebulizer and cyclonic spray chamber, and a 10-second integration.

Element	Wavelength (nm)	D.L.
Al	396.152	0.9
As	193.696	3.6
B	249.772	0.25
Be	313.107	0.017
Cd	214.440	0.07
Co	238.892	0.25
Cr	267.716	0.25
Cu	224.700	0.9
Fe	259.939	0.2
Mn	257.610	0.03
Mo	202.031	2
Ni	231.604	0.4
Pb	220.353	1.4
Sb	206.836	4
Se	196.026	4.5
Tl	190.801	3.5
V	309.310	0.15
Zn	206.200	0.2

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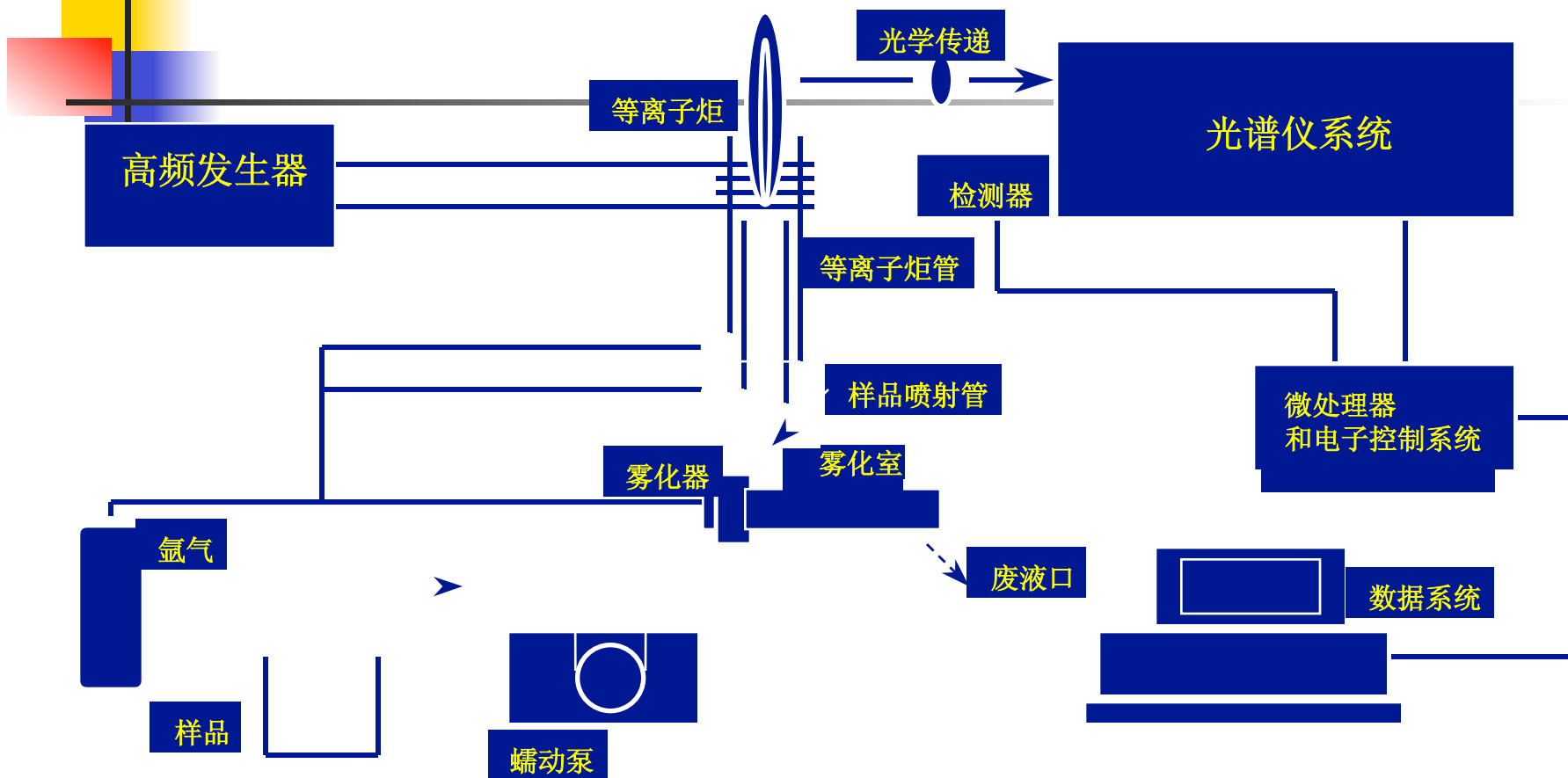
Suite 5401, Central Plaza, No.18 Harbour Road, HONG KONG

Optima ICP-OES, Advantage and Disadvantage

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Suite 5401, Central Plaza, No.18 Harbour Road, HONG KONG

AAnalyst 400 with MHS-15

- Low cost comparing to Optima ICP-OES
- Easy to use
 - WinLab32 software
 - Computer-controlled gas box (C₂H₂ + Compressed Air)
- Easy to maintain and service
 - Quick-lock Burner Assembly
 - User-replaceable parts...
- Performance
 - Daily sample throughput 25-30 samples

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Other similar techniques:

得利環球有限公司 Atomic Absorption Spectroscopy (AAS) Tel: 2802-3322 Fax: 2802 8988

Inductively Coupled Plasma Optical Emission Spectroscopy (ICP-OES)

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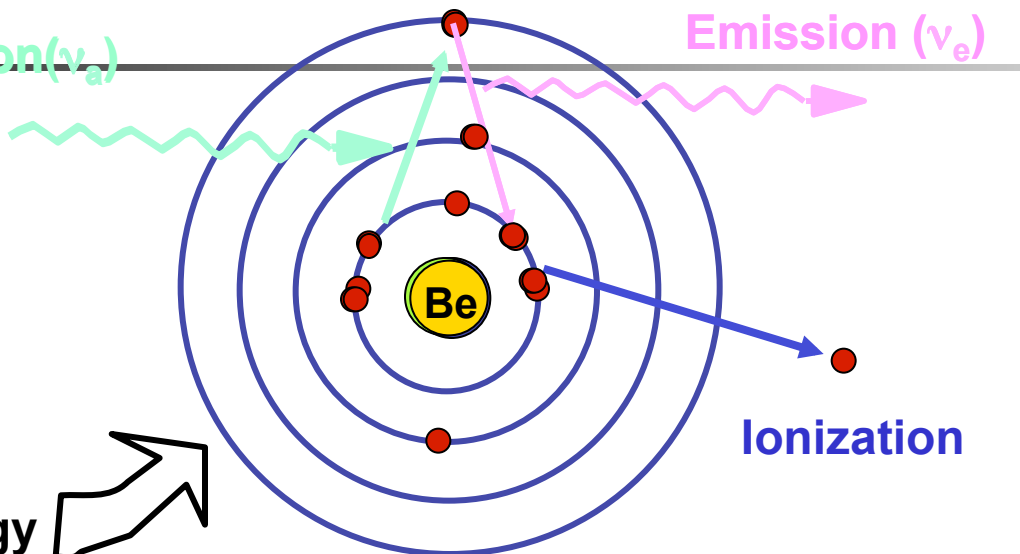
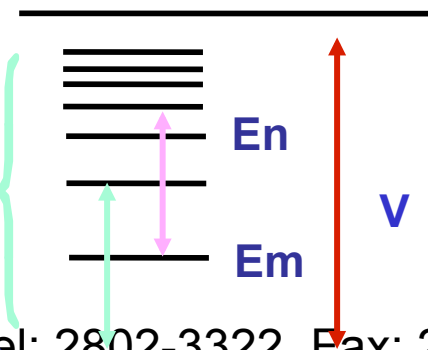
Suite 5401, Central Plaza, No.18 Harbour Road, HONG KONG

ICP-MS 技術

Advantages

- Exceptional detection limits
10⁻¹² 数量级
- Multiple element technique
多元素分析
- Excellent dynamic range
动态范围宽
- High productivity
速度快
- Integration of alternate sampling

Limitations

- Expensive to purchase
价格贵
- Training time to gain skill
仪器操作相对复杂

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- The Optima 2100DV
 - Typically less than 100 samples a day (每天分析100个样品)
 - 1 to 10 elements per sample (10个元素/样品)
 - Best analytical performance
- The Optima 5300DV
 - Challenging analytical problems (多元素/样品, 复杂基体)
 - Demands sample throughput (每天分析150个以上样品)
 - Ultimate analytical performance
- The AAnalyst 400
 - 30 samples per day (每天分析大约30个样品)

Installation Requirement: Lab Space and Bench

Heavy duty (over 200kg) and anti-corrosive bench for Optima ICP-OES & AAS

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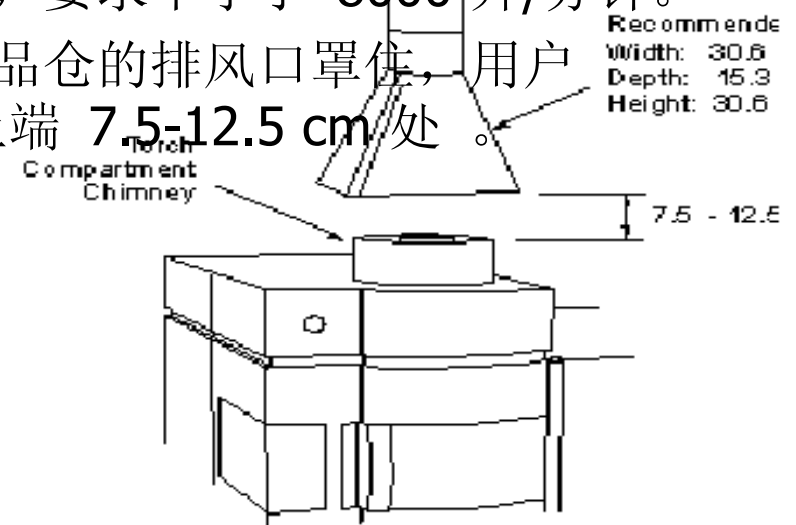
Photo: The Hong Kong University of Science and Technology

Installation Requirement: Ventilation

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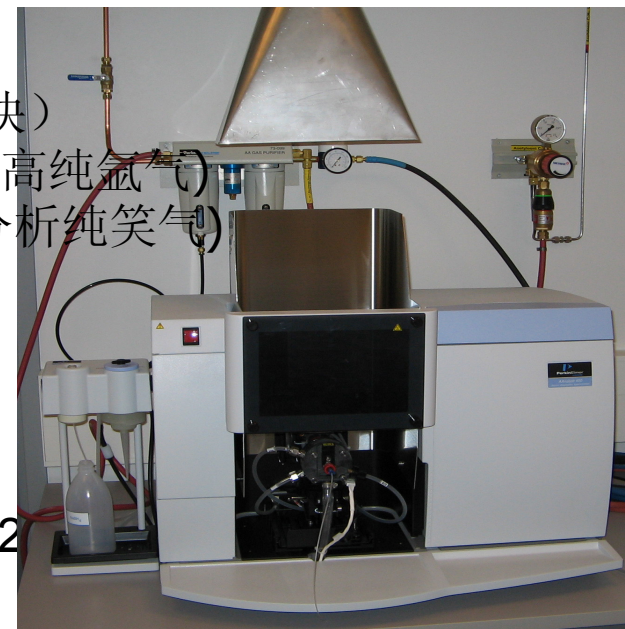
Installation Requirement: Electricity

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Installation Requirement: Gas (高纯氩气)

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- VAT (资本增值税, in RMB) : 17% Contract Price,
- optional: 2% handling charge for import fees, custom clearance and etc.
- Tariff (海关关税): ZERO
- Custom Code: 9027 3000
- IPPC (木质包装认证) : provided by PerkinElmer (Hong Kong) Limited

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